



Maleimide-Activated Ovalbumin

Applicable to:

407-0002 2 mg lyophilized maleimide-activated ovalbumin

407-0005 5 mg lyophilized maleimide-activated ovalbumin

Release4

29/06/2016

Introduction

Maleimide-activated Ovalbumin is designed to facilitate the conjugation of ovalbumin to proteins, peptides and ligands that contain sulfhydryl (-SH) groups. Ovalbumin stimulates a large immune response and is typically used as a carrier protein to generate antibodies against small antigens (peptides or organic compounds). For success the antigen (normally referred to as a hapten) needs to be covalently linked to the carrier protein. The added maleimide groups allow conjugation to be performed very efficiently at physiological pH.

Storage and shipping

Maleimide-activated ovalbumin is shipped at ambient temperature as a lyophilized powder and should be stored upon receipt at -20°C. Shelf life is 3 years.

Technical notes

The buffer in which the thiolated molecule is dissolved should be in the pH range 6.5-7.5. Phosphate buffered saline pH 7.2 is commonly employed. EDTA (1-10 mM) may help to prevent metal catalysed oxidation of thiols. Please see the below table for recommended buffer conditions and components:

Buffer components	
pH range 6.5-7.5	✓
Chelating agents (e.g. EDTA)	✓
Primary amines	✓
Reducing reagents	X
Sodium Azide	< 0.1%

Molecules that do not contain -SH groups can be chemically modified to introduce the required thiol functionality. Innova offers a thiolation kit (Prod. Code: 419-0005) compatible with this product. In the case of antibodies, disulfide (S-S) bridges may also be cleaved to release free thiols. Any reducing reagent (e. g. dithiothreitol and β -mercaptoethanol) used to produce free thiols should be removed before proceeding with the conjugation. Further technical information about chemical modification strategies is available on our website.

Limited use license

Innova Biosciences' Maleimide-activated labels are offered for research purposes alone, and are not intended for human, therapeutic or diagnostic use. The purchase of this product conveys to the buyer (whether the buyer is a not-for-profit, academic or for-profit entity) the non-transferable right to use the amount of product purchased and the components of the product for in-house research. The buyer shall not sell or otherwise transfer this product, its components, or materials prepared therefrom to any third party. The buyer shall not use this product or its components for commercial purposes. For the avoidance of doubt, 'commercial purposes' means any activity by a party for consideration and includes, without limitation, use of the product or its components (i) in the manufacturing of conjugated materials (e.g. labeled antibodies), (ii) to provide a service, information or data, (iii) for therapeutic, diagnostic or prophylactic purposes, or (iv) for repackaging/resale, whether or not such product or its components are resold for use in research. The use of this product by the buyer constitutes agreement with the terms of this limited use label license for Innova Biosciences' products. For information on purchasing a license for commercial applications contact Innova Biosciences Ltd, Business Development Office, Babraham Hall, Babraham, Cambridge, UK, CB22 3AT. Tel +44(0)1223 661000; Fax +44(0)1223 496172.

Protocol

1. Allow the vial of maleimide-activated ovalbumin to warm to room temperature.
2. Remove the crimp seal and reconstitute the lyophilized powder by adding a buffered solution of the thiol-containing molecule. Gently agitate the vial to ensure thorough dissolution.
3. Incubate for 3-4 hours (or overnight) at room temperature and then store at 4°C.

Related products

Description	Prod. Code
Maleimide Alkaline Phosphatase	402-0005
Maleimide HRP	401-0005
Maleimide Streptavidin	405-0005
Maleimide R-Phycoerythrin	403-0005
Maleimide APC	404-0002
Maleimide BSA	408-0005
Maleimide KLH	409-0005
Thiolation kit	419-0005
Thiol quantification kit	418-0002

For further information about our products please see our website: www.innovabiosciences.com.

For technical support please contact us: www.innovabiosciences.com/contact-us.html